

Cyto System

Order No.:
Cyto System

Rotor:
C0326-72

Buckets:
C0323-010

Cyto container:
C0323-064
cylindrical, 2 x 1.5mL
C0323-065
conical, 2 x 3mL

Carrier:
C0323-062

Filter cards:
C0323-085 (cylindrical)
C0323-086 (conical)

Lid:
C0323-103

Sedimentations occur at rcf-values
between 40 and 690 x g
This is equivalent to speeds between
530 and 2110 rpm

Carrier



Filter card



Cyto container



Micro tubes



Buckets



Rotor



Lid



1 Step

Insert the microscope slide
into the carrier

2 Step

Place a filter card onto the
microscope slide

3 Step

Put the cyto container onto
the filter card.

4 Step

Place the micro tube (with
bottom hole) into the
container chamber, and fill it
with the sample fluid

5 Step

Insert the carriers including the
container into the bucket

6 Step

Put the buckets into the
centrifuge rotor immediately

7 Step

If necessary, close the
container with the lid

Application possibilities

Z 306

Z 326

Z 326K

Z 306	310.00 V01 - Z 306, 230V / 50-60 Hz / 240W 310.00 V02 - Z 306, 120V / 50-60 Hz / 300W
Z 326	312.00 V01 - Z 326, 230V / 50-60 Hz / 240W 312.00 V02 - Z 326, 120V / 50-60 Hz / 300W
Z 326K	311.00 V01 - Z 326K, 230V / 50-60 Hz / 240W 311.00 V02 - Z 326K, 120V / 50-60 Hz / 300W

